Protein G Terbium Fluorescence Magnetic Beads

BcMagTM Protein G Terbium Fluorescence Magnetic Beads (Fig.1) are Time-Resolved Fluorescence (TRF) magnetic microspheres covalently coupled with Protein G on the surface. The beads are manufactured using nanometer-scale superparamagnetic iron oxide and terbium metal as core and entirely encapsulated by a high-purity silica shell, ensuring no leaching problems with the iron oxide and europium metal. The microspheres combine the benefits of a novel avidin biotin-binding method, time-resolved Fluorescence dyes, and magnetic characteristics to perform very sensitive assays.



Fig.1 BcMagTM Terbium Fluorescence Magnetic Beads

The recombinant Protein G contains 200 amino acids having a molecular mass of 21.8kDa with two Fc-binding domains per protein. These domains can bind to antibodies from many species, including mice, humans, rabbits, cows, goats, and sheep. The albumin and cell surface binding domains have been removed from Recombinant Protein G to reduce nonspecific binding. This recombinant protein has a significantly greater capacity and more specific IgG binding than native Protein G and Protein A for most mammalian IgGs. It can purify mammalian IgGs that do not bind well to Protein A. In addition, when working with complicated biological samples, the beads blocked by our unique blocking agent minimize and reduce nonspecific binding to the bead surface. Table 1 lists the binding properties.

Species	Antibody	Binding	Species	Antibody	Binding
		(Protein G)			(Protein G)
	lgG 1	+++		lgG1	++++
	IgG 3	++++	Sheep	IgG2	++++
Mouse	IgG 2a	++++		Total IgG	++++
	IgG 2b	++++		IgG(ab)	-
	IgM	-		IgG(c)	-
	Total IgG	++++	Horse	IgG(T)	++++
	lgG1	++++		Total IgG	++++
	IgG2	++++		lgG1	++++
	IgG3	++++	Goat	IgG2	++++
	IgG4	++++		Total IgG	++++
Human	IgA	-		lgG1	++++
	IgD	-	Cow	IgG2	++++
	IgM	-		Total IgG	++++
	Fab	++	Rabbit	Total IgG	++++
	scFv	-	Guinea Pig	Total IgG	++
	Total IgG	++++	Pig	Total IgG	++
	lgG 1	+++	Cat	Total IgG	++
	IgG 2a	++++	Dog	Total IgG	++



Magnetic Beads Make Things Simple

Instruction Manual

2

Rat	IgG 2b	++	++++ (Strong Binding); +++ (Medium Binding);
	IgG 2c	++++	++ (Weak Binding); - (No Binding); N/A (No
	Total IgG	+++	Information)
			Table 1. Protein G binding properties

Although conventional fluorophores have been widely used over the past decades, they still suffer from either one or several limitations in terms of applicability and efficiency: 1. Narrow excitation bands cause higher background signals. 2. Smaller Stokes shift often produces self-quenching. 3. Fluorescence is sensitive to environmental factors such as metallic ion concentration, pH, temperature, and solvent polarity. 4. Fluorescence intensity is not high enough for detecting a single biomolecular. 5. Fluorescence intermittency (blinking) affects some processes of molecule detection. 6. Easily aggregated because of hydrophobicity.

Fluorescence dye properties

	Fluorophore	Fluorescence color	Excitation (nm)	Emission (nm)	Fluorescence lifetime (T) (µsec)	Stokes shifts (nm)	Selection of Emission Filter
Ī	Terbium (Tb ³⁺)	Green	320	545	1050	220	545/40

Specification			
Bead Size	2.5μm diameter; 5μm diameter		
Number of Beads	~10 x 10 ⁷ beads/mg (2.5μm), ~5 x 10 ⁷ beads /mg (5μm)		
Magnetization	~40-45 EMU/g		
Type of Magnetization	Superparamagnetic		
Concentration	10 mg/ml (10mM Tris, 0.15 M NaCl,0.1% BSA,1 mM EDTA,		
	pH7.4)		
Binding Capacity	~ 1 mg lgG/ml of Beads		
Storage	Ship at room temperature. Store at 4°C. Do not freeze		

BcMagTM TR-FRET (Time-Resolved FRET) Assay

BcMagTM TR-FRET Assay, in contrast to typical FRET (Förster Resonance Energy Transfer) assays, uses time-resolved Fluorescence magnetic beads (BcMagTM TR-Magnetic Beads) as the donor fluorophore. The donor and acceptor can be two proteins, two DNA strands, an antigen, an antibody, or a ligand and its receptor. After a reasonable time delay (usually 50 to 100 s), a signal is generated by fluorescence resonance energy transfer between a donor and an acceptor molecule when they are close and monitored in a time-resolved way. In BcMagTM TR-FRET Assay, a trace amount of analytes can be easily enriched from the complex by TR-Magnetic Beads, resulting in higher sensitivity. This assay practically eliminates all fluorescence backgrounds caused by the sample and plastic microplate, as well as by direct acceptor excitation. As a result, the signal-to-noise ratios of the BcMagTM TR-FRET Assay are very high, and the background is quite low. Furthermore, the assay does not need washing steps.

BcMagTM TR-FRET Assay offers substantial advantages to bioassays in high throughput screening, such as assay flexibility, dependability, increased assay sensitivity, higher throughput, and fewer false positive/false negative results.

Terbium cryptate fluorophore is an efficient Fluorescence label due to its distinct specific properties. It is excited at 320nm and emits green fluorescence at 545nm, with a long fluorescence lifetime (1050 μsec) and large stokes shifts (220 nm). By taking advantage of these properties, time-resolved fluorescence measurement can dramatically reduce the fluorescence background from the sample and increase the signal-to-noise ratio to offer detectability better than one order of magnitude than conventional Fluorescence dyes. BcMagTM Protein A/G Terbium Fluorescence Magnetic Beads are excellent donors used in TR-FRET assays.

Workflow of TR-FRET Magnetic Beads Assay (Fig.2)

1. Mix the antibody-conjugated donor beads with the cell lysates and incubate them with continuous rotation for a sufficient time. The beads remain suspended in the sample solution during mixing, allowing the target analytes to bind to the donor beads. 2. After incubation, the beads are collected and separated from the sample using a magnet rack. 3. Add the antibody-conjugated acceptor and incubate them with continuous rotation for a sufficient time. 4. Analysis of numerous microplate readers supports TR-FRET measurements.



Magnetic Beads Make Things Simple

Instruction Manual

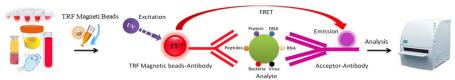


Fig.2 Workflow of TR-FRET Magnetic Beads Assay

Advantages and benefits

- 1. Perform a double function simultaneously on the same beads: The magnetic beads combine separation/preconcentration and detect analytes, allowing quick, simple, robust, and high-throughput analytes of trace amounts from complex biological samples on the same beads.
- 2. Ultra sensitive. Lower detection limits of 10 pg/mL versus typical fluorometric detection limits of 100 pg/mL
- 3. Extremely photostable and highly resistant to photobleaching. All the lanthanide chelate or cryptate molecules and iron oxide are entirely encapsulated inside each bead instead of merely on the bead's surface. The protective environment prevents iron oxide and dye from leaching into aqueous media, which makes the beads less sensitive to external conditions such as solvent, temperature, pH, etc.
- 4. Very high Fluorescence intensity. Because a single bead has a large concentration of lanthanide chelate with a high quantum yield ranging from 40 to 90%, the beads show excellent fluorescence intensity, which increases test sensitivity without signal amplification. Such bright beads are also perfect for donors' use in time-resolved FRET assays.
- Lanthanide chelate or cryptate has large Stokes shifts (>250 nm), narrow emission bands (-10 nm bandwidth), and long fluorescence lifetime
 (μs), which dramatically reduces background and increases the signal-to-noise ratio.
- 6. Most bioprocess ELISA assays can be converted to an HTRF assay.
- 7. No washing step is involved in the assays.
- 8. Have a hydrophilic silica surface grafted by different functional groups with linkers of variable lengths, allowing efficient conjugation of various ligands such as peptides, proteins, antibodies, small molecules, carbohydrates, aptamers, DNA/RNA, etc.
- 9. Due to the microsphere's magnetic properties, the Fluorescence magnetic beads are suitable for high-throughput automation.

Related Products				
Streptavidin Europium Fluorescent Magnetic Beads	Aldehyde-Activated Europium Fluorescent Magnetic Beads			
Streptavidin Terbium Fluorescent Magnetic Beads	Aldehyde-Activated Terbium Fluorescent Magnetic Beads			
Streptavidin-Ruthenium Fluorescent Magnetic Beads	Aldehyde-Activated Ruthenium Fluorescent Magnetic Beads			
Avidin Europium Fluorescent Magnetic Beads	Amine Activated-Europium Fluorescent Magnetic Beads			
Avidin Terbium Fluorescent Magnetic Beads	Amine-Activated Terbium Fluorescent Magnetic Beads			
Avidin Ruthenium Fluorescent Magnetic Beads	Amine-Activated Ruthenium Fluorescent Magnetic Beads			
Protein A and G Europium Fluorescent Magnetic Beads	Carboxyl-Activated Europium Fluorescent Magnetic Beads			
Protein A and G Terbium Fluorescent Magnetic Beads	Carboxyl-Activated Terbium Fluorescent Magnetic Beads			
Protein A and G Ruthenium Fluorescent Magnetic Beads	Carboxyl-Activated Ruthenium Fluorescent Magnetic Beads			
Protein A Europium Fluorescent Magnetic Beads	Hydrazide-Activated Europium Fluorescent Magnetic Beads			
Protein A Terbium Fluorescent Magnetic Beads	Hydrazide-Activated Terbium Fluorescent Magnetic Beads			
Protein A Ruthenium Fluorescent Magnetic Beads	Hydrazide-Activated Ruthenium Fluorescent Magnetic Beads			
Protein G Europium Fluorescent Magnetic Beads	Iodoacetyl-Activated Europium Fluorescent Magnetic Beads			
Protein G Terbium Fluorescent Magnetic Beads	Iodoacetyl-Activated Terbium Fluorescent Magnetic Beads			
Protein G Ruthenium Fluorescent Magnetic Beads	Iodoacetyl-Activated Ruthenium Fluorescent Magnetic Beads			
Protein L Europium Fluorescent Magnetic Beads	NHS-Activated Europium Fluorescent Magnetic Beads			
Protein L Terbium Fluorescent Magnetic Beads	NHS-Activated Terbium Fluorescent Magnetic Beads			



Magnetic Beads Make Things Simple

Instruction Manual

Protein L-Ruthenium Fluorescent Magnetic Beads	NHS-Activated Ruthenium Fluorescent Magnetic Beads
--	--